

SSR-3A-MC Autoclave Quick Guide

Please see procedure Autoclave SSR-3A-MC Use and Operation [EQUIP 2.0] for full instructions, safety information, error messages, and troubleshooting.

1. Fill out Autoclave Use Log (on fridge across from autoclave).
2. If screen is blank, press any key to turn it on.
3. Wearing proper PPE, load autoclave.
 - a. Ensure all items are in buckets and all caps/closures are loose.
4. Press Jacket button. "JACKET PRESSURE IS ON" will display on top right corner of screen.
5. From the MAIN MENU, select the cycle (FAST, DRY or FLUID) using corresponding keypad number.
 - a. If MAIN MENU is not displayed, press 9 to exit until MAIN MENU screen appears.
6. Press 9 twice to return to the MAIN MENU.
7. Press RUN button on keypad.
8. Press ENTER twice to begin run.
 - a. It is not necessary to enter load or user number.
9. After run is complete, **make sure jacket pressure is set to OFF.**
 - a. Jacket will turn itself back on after DRY cycle!